



LabStrong Cascade DI Type II Water System

Cartridge Placement Order

Dual Style Cartridge Systems (D00292)

Instructions:

1. Remove cartridges from plastic bag and remove sealing plugs at each end of the cartridge
2. Place cartridges in Cascade holder in the order identified in the chart below and as shown in figure 1
3. Tighten Velcro straps around cartridges
4. Once all cartridges are in place, ensure tubing is securely locked into position as shown in figure 2.
5. Securely connect fittings at top and bottom of each cartridge and attach inlet water to system .
6. Flow rate should be 1.0-1.5 Liters per minute with a max of 1.8 Liters per minute.

Please refer to complete Cascade manual (LT 0029-1) provided with the LabStrong Cascade Water System for detailed information. **Scan the QR code for how-to video of installation!**

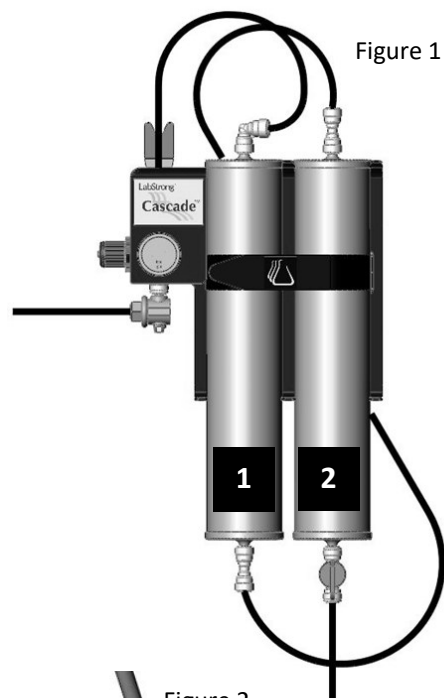


Figure 1

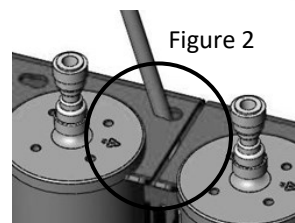


Figure 2



Dual Style Cartridge Systems

D002912KT Pretreat Organic		D002913KT Pretreat Standard		D002914KT Organic Removal (carbon)	
Part #	Description	Part #	Description	Part #	Description
D00291	PRETREAT ORGANIC / CARBON	D00291	PRETREAT ORGANIC / CARBON	D00298	ORGANIC REMOVAL (CARBON)
D00296	ORGANIC FREE (Dyed Resin)	D00297	ULTRAPURE (Dyed Resin)	D00298	ORGANIC REMOVAL (CARBON)

D002916KT Pretreat Distillation	
Part #	Description
D00291	PRETREAT ORGANIC / CARBON
D002915	HIGH CAPACITY (DYED RESIN)

LabStrong Corporation
7709 Commerce Park
Dubuque, Iowa 52002

Phone: (563) 588-8900
Fax: (563) 588-3601
Sales@labstrong.com

www.LabStrong.com





LabStrong Cascade DI Type II Water System


Cartridge Placement

Quad Style Cartridge Systems (D00294)



Instructions

1. Remove cartridges from plastic bag and remove sealing plugs at each end of the cartridge
2. Begin placing cartridges in Cascade holder in the order identified in the chart below and as shown in Figure 1
3. Tighten Velcro straps around cartridges
4. Once all cartridges are in place, ensure tubing is securely locked into position as shown in Figure 2.
5. Securely connect fittings at top and bottom of each cartridge and attach inlet water to system.
6. Flow rate should be 1.0-1.5 Liters per minute with a max of 1.8 Liters per minute.

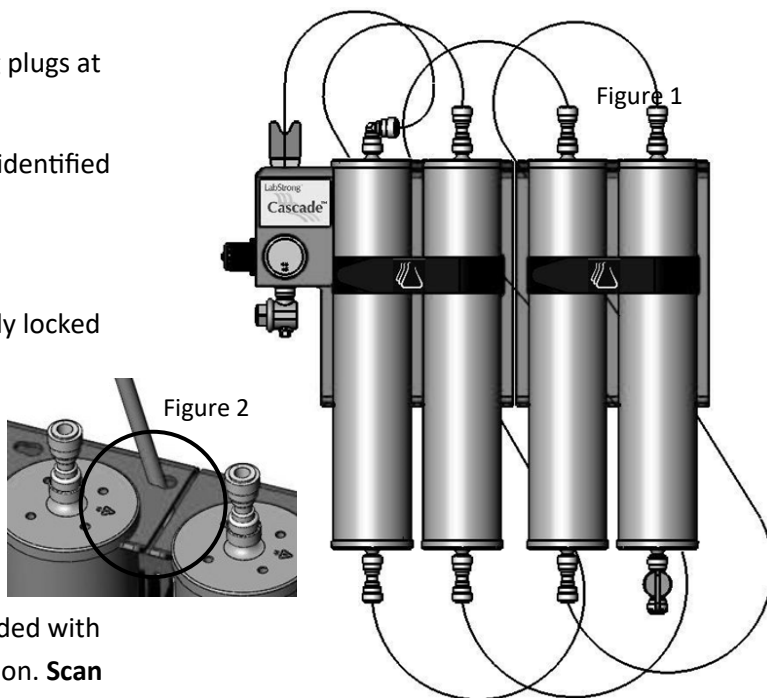

Please refer to complete Cascade manual (LT 0029-1) provided with the LabStrong Cascade Water System for detailed information. **Scan the QR code for how-to video of installation!**

Quad Style Cartridge Systems

Tap Water Feed

D00298KT Organic Kit		D00299KT Standard Kit	
Part #	Description	Part #	Description
1 D00291	PRETREAT ORGANIC/CARBON	D00291	PRETREAT ORGANIC/CARBON
2 D00293	HIGH CAPACITY	D00293	HIGH CAPACITY
3 D00295	ULTRAPURE	D00295	ULTRAPURE
4 D00296	ORGANIC FREE (Dyed Resin)	D00297	ULTRAPURE (Dyed Resin)

Pretreat Water Feed

D002910KT Organic Pretreat Kit		D002911KT Standard Pretreat Kit	
Part #	Description	Part #	Description
1 D00291	PRETREAT ORGANIC/CARBON	D00291	PRETREAT ORGANIC/CARBON
2 D00295	ULTRAPURE	D00295	ULTRAPURE
3 D00295	ULTRAPURE	D00295	ULTRAPURE
4 D00296	ORGANIC FREE (Dyed Resin)	D00297	ULTRAPURE (Dyed Resin)